Safety Information for Genomics Instruments

For use with Access Array™, Biomark™/Biomark HD, C1™, Callisto™, EP1™, FC1™, IFC Controllers, Juno™, and Polaris™

Contents

Safety Guide	2
Introduction	2
Safety Guidelines	3
Sikkerhedsvejledning	7
Indledning	7
Sikkerhedsretningslinjer	8
Sicherheitsleitfaden	12
Einleitung	12
Sicherheitsleitlinien	13
Guía de seguridad	17
Introducción	17
Pautas de seguridad	18
Guide de sécurité	22
Introduction	22
Directives de sécurité	23
Guida alla sicurezza	27
Introduzione	27
Linee guida sulla sicurezza	28
Veiligheidshandleiding	32
Inleiding	32
Veiligheidsvoorschriften	33

Safety Guide

For Research Use Only. Not for use in diagnostic procedures.

Introduction

This document provides important safety information pertaining to the operation of Fluidigm genomics instruments. For detailed information about using an instrument, see the appropriate user guide.

IMPORTANT Before using the instrument, read and understand the safety guidelines in this document. Failure to follow these guidelines may result in undesirable effects, injury to personnel, and/or damage to the instrument or to property.

Safety Alert Conventions

Fluidigm documentation uses specific conventions for presenting information that may require your attention. Refer to the following safety alert conventions.

Safety Alerts for Chemicals

For hazards associated with chemicals, this document follows the United Nations Globally Harmonized System of Classification and Labelling of Chemicals (GHS) and uses indicators that include a pictogram and a signal word that indicates the severity level:

Indicator	Description
	Pictogram (see example) consisting of a symbol on a white background within a red diamond-shaped frame. Refer to the individual safety data sheet (SDS) for the applicable pictograms and hazards pertaining to the chemicals being used.
DANGER	Signal word that indicates more severe hazards.
WARNING	Signal word that indicates less severe hazards.

Safety Alerts for Instruments

For hazards associated with instruments, this document uses indicators that include a pictogram and signal words that indicate the severity level:

Indicator	Description
	Pictogram (see example) consisting of a symbol on a white background within a black triangle-shaped frame. Refer to the instrument user guide for the applicable pictograms and hazards pertaining to instrument usage.
DANGER	Signal word that indicates an imminent hazard that will result in severe injury or death if not avoided.
WARNING	Signal word that indicates a potentially hazardous situation that could result in serious injury or death if not avoided.
CAUTION	Signal word that indicates a potentially hazardous situation that could result in minor or moderate personal injury if not avoided.
IMPORTANT	Signal word that indicates information necessary for proper use of products or successful outcome of experiments.

Safety Data Sheets

Read and understand the SDSs before handling chemicals. To obtain SDSs for chemicals ordered from Fluidigm, either alone or as part of this system, go to fluidigm.com/sds and search for the SDS using either the product name or the part number.

Some chemicals referred to in this user guide may not have been provided with your system. Obtain the SDSs for chemicals provided by other manufacturers from those manufacturers.

Safety Guidelines

General Safety

In addition to your site-specific safety requirements, Fluidigm recommends the following general safety guidelines in all laboratory and manufacturing areas:

- Use the appropriate personal protective equipment (PPE): safety glasses, fully enclosed shoes, lab coats, and gloves, according to your laboratory safety practices.
- Know the locations of all safety equipment (fire extinguishers, spill kits, eyewashes/showers, first-aid kits, safety data sheets, etc.), emergency exit locations, and emergency/injury reporting procedures.
- Do not eat, drink, or smoke in lab areas.
- Maintain clean work areas.
- Wash hands before leaving the lab.

Instrument Safety

The instrument should be serviced by authorized personnel only.

WARNING Do not modify this instrument. Unauthorized modifications may create a safety hazard.

WARNING BIOHAZARD. If you are putting biohazardous material on the instrument, use appropriate personal protective equipment and adhere to Biosafety in Microbiological and Biomedical Laboratories (BMBL), a publication from the Centers for Disease Control and Prevention, and to your lab's safety protocol to limit biohazard risks. If biohazardous materials are used, properly label the equipment as a biohazard. For more information, see the BMBL guidelines online at cdc.gov/biosafety/publications/index.htm.

WARNING PHYSICAL INJURY HAZARD. Do not attempt to lift or move any boxed or crated items unless you use proper lifting techniques. See the user guide for the weight of the boxed or crated instrument.

If you choose to lift or move the instrument after it has been installed, do not attempt to do so without the assistance of others. Use appropriate moving equipment and proper lifting techniques to minimize the chance of physical injury.

CAUTION PINCH HAZARD. The instrument door and tray can pinch your hand. Make sure your fingers, hands, and shirtsleeves are clear of the door and tray when loading or ejecting an integrated fluidic circuit (IFC).

CAUTION HOT SURFACE HAZARD. The thermal cycler chuck gets hot and can burn your skin. Use caution when working near the chuck.

CAUTION HOT SURFACE HAZARD. Never press down on the integrated fluidic circuit (IFC) when it is on the thermal cycler chuck. If you encounter a vacuum problem, turn off the system, allow it to cool down, and remove the IFC. Clean the bottom of the IFC and/or chuck surface with a lint-free cloth and 70% isopropyl alcohol.

CAUTION HOT SURFACE HAZARD. Make sure the chuck has had time to cool. It can get very hot and cause burn injury.

CAUTION HOT SURFACE HAZARD. Be careful not to touch the heated anvil on the IFC Controller RX and IFC Controller WX after a purge script has been run or is aborted. The anvil is normally not hot enough to cause injury, but if the integrated fluidic circuit (IFC) is ejected during a purge script, the anvil may be as hot as 90 °C.

Symbols on the Instrument

The following table describes the hazard symbols that may be used in this document or on labels on the instrument.

Symbol	Description
	Hazard. Consult the user guide for further information.
	Hot surface hazard. Do not touch; potential for personal injury.
	Biohazard.
	Electricity hazard. Indicates high electricity levels and a threat of electric shock from machines and/or equipment in the vicinity. You may suffer severe injuries or death.
	Pinch hazard. Indicates where pinch hazards exist. Exercise caution when operating around these areas.
	
	
	Lifting hazard.
	Indicates specific chemical harm.
	Indicates hazardous, toxic, or very toxic materials that are very hazardous to health or potentially fatal when inhaled, swallowed, or in contact with the skin.
	Indicates a health hazard.
	Power and standby symbol.
	Power switch is in the Off position.
	Power switch is in the On position.
	Protective conductor terminal (main ground). It must be connected to earth ground before any other electrical connections are made to the instrument.
	To minimize negative environmental impact from disposal of electronic waste, do not dispose of electronic waste in unsorted municipal waste.
	Follow local municipal waste ordinances for proper disposal provision. Contact customer service for information about responsible disposal options.

Electrical Safety

NOTE The main power disconnect is on the rear panel of the instrument.

WARNING ELECTRICAL HAZARD. DO NOT REMOVE THE COVERS. Electrical shock can result if the instrument is operated without its protective covers. No internal components are serviceable by the user.

WARNING ELECTRICAL HAZARD. Plug the instrument into a properly grounded receptacle with adequate current capacity.

Chemical Safety

The responsible individuals must take the necessary precautions to ensure that the surrounding workplace is safe and that instrument operators are not exposed to hazardous levels of toxic substances. When working with any chemicals, refer to the applicable safety data sheets (SDSs) provided by the manufacturer or supplier.

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For technical support visit fluidigm.com/support.

North America +1 650 266 6100 | Toll-free (US/CAN): 866 358 4354 | techsupport@fluidigm.com Latin America +1 650 266 6100 | techsupportlatam@fluidigm.com
Europe/Middle East/Africa/Russia +44 1223 859941 | techsupporteurope@fluidigm.com China (excluding Hong Kong) +86 21 3255 8368 | techsupportchina@fluidigm.com
Japan +81 3 3662 2150 | techsupportjapan@fluidigm.com All other Asian countries/India/Australia +1 650 266 6100 | techsupportasia@fluidigm.com

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Information in this publication is subject to change without notice. Patent and license information: fluidigm.com/legalnotices. Fluidigm, the Fluidigm logo, Access Array, Biomark, C1, Callisto, EP1, FC1, Juno, and Polaris are trademarks or registered trademarks of Fluidigm Corporation in the United States and/or other countries. © 2018 Fluidigm Corporation. All rights reserved. 04/2018